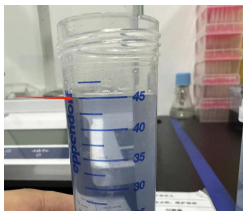
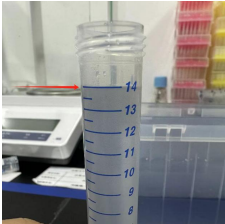


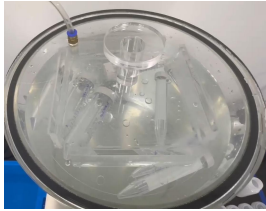

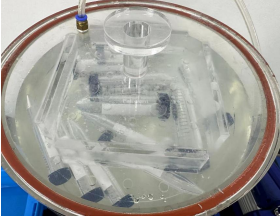







Eppendorf & PakGent Centrifuge tubes compare report

Product:	15/50ML centrifuge tubes	quantity:	6 pcs each model	test equipment:	centrifugal machine seal tester autoclave ultra low temperature refrigerator	test date:	2024/7/9-7/10
test item	requirement	test photo				conclusion	
		Eppendorf		PakGent		Eppendorf	PakGent
calibration test	allow +/-5% tolerance					Pass the test	Pass the test
seal test	70kpa, 20 seconds ,no leakage		 艾本德离心管密封测试.mp4		 莱杰离心管密封测试.mp4	15/50ml centrifuge tube pass the test	15/50ml centrifuge tube pass the test
autoclave test	First autoclave the 50 CT tubes and cool it Add 80% water, 15000g, 10 min, no broken and leakage					<ul style="list-style-type: none"> 1.15/50ml centrifuge tube no sharp change 2.15/50ml centrifuge tube no liquid leakage 3.15/50ml centrifuge tube writing area no tear 4.15/50ml centrifuge tubes have small cracks on the bottom 	<ul style="list-style-type: none"> 1.15/50ml centrifuge tube no sharp change 2.15/50ml centrifuge tube no liquid leakage 3.15/50ml centrifuge tube writing area no tear 4.15/50ml centrifuge tubes no cracks on the tube body and bottom

<p>cryopreserved test</p>	<p>First put the CT tubes at -80°C, 24hrs, after the tubes warm to room temp Add 80% water, 15000g, 10 min, no broken and leakage</p>	   	   		<p>1.15/50ml centrifuge tube no sharp change 2.15/50ml centrifuge tube no liquid leakage 3.15/50ml centrifuge tube writing area no tear 4.15/50ml centrifuge tubes have small cracks on the bottom</p>	<p>1.15/50ml centrifuge tube no sharp change 2.15/50ml centrifuge tube no liquid leakage 3.15/50ml centrifuge tube writing area no tear 4.15/50ml centrifuge tubes no cracks on the tube body and bottom</p>
---------------------------	---	---	---	---	---	--